

Wash Baskets & Machine Accessories

For use in Miele Medium Chamber Washers Models G 7825 & G 7826



Miele Baskets & Accessories: Advancing the Science of Clean

Engineered Simplicity

Miele washers expertly clean a wide variety of items through a simple-to-use system of interchangeable baskets and inserts. Miele wash baskets are engineered to roll directly onto the transfer cart, with a modular basket design allowing simple removal of each level away from the machine. This ergonomic design allows easy loading and unloading at the washer or at the glassware storage cabinet.

This interchangeable basket system ensures a precise match to your specific cleaning requirements. And, thanks to Miele's universal design for medium chamber wash baskets and inserts, a new basket will fit a new Miele G 7825 washer as precisely as it will an older model.

To further simplify basket and wash cycle selection, the G 7825 features an optional Automatic Basket Recognition, ensuring repeatable wash results and eliminating guesswork from the machine operator regarding wash parameters.

Injection Cleaning and Drying

The Miele G 7825 washer features 1 to 3 levels of injection cleaning or up to 5 levels of non-injection cleaning, providing excellent throughput and wide array of wash load possibilities.

Miele wash baskets feature removable and interchangeable injectors for a truly custom basket configuration. This interchangeability also allows easy replacement of older injectors, ensuring many years of useful operation.

When Miele's HEPA-filtered forced air drying is selected, air is circulated throughout the wash chamber and through the injectors to the interior of narrow-neck flasks. This provides superior drying typically within 15 to 30 minutes, eliminating the need for a separate drying oven. This system frees up valuable floor space, reduces labor costs, and lowers your electrical costs.

Gentle Yet Effective Cleaning

Miele washers utilize high-volume, low-pressure water circulation to prevent breakage inside the wash chamber.

This care is carried over into our wash basket design. Each basket delicately cradles the item being washed while providing renowned, repeatable wash results. An excellent example is pipettes. Miele provides perfect results for difficult items like this, eliminating breakage even for small items.

Solutions for Your Specific Application

If you can not find a basket solution, please contact us for more information. Through the exclusive Miele Application Laboratory, housed in Princeton, NJ, with a sister lab in Gütersloh, Germany, we are continually developing new wash baskets, ensuring you a competent, customized solution to your cleaning requirements.

Mobile Baskets



E 775 Mobile Unit

- With 2 levels for various inserts
- Built-in middle spray arm
- Magnetic coding strip for automatic mobile unit sensor (optional)
- Includes 2 E 750 units to secure and position inserts for optimal cleaning results
- Clearance: (Level 1 = Lower Level)
1: 304 mm H, 482 mm W, 590 mm D
2: 290 mm H, 488 mm W, 546 mm D
- Dimensions: Height 400 mm
Width 530 mm
Depth 600 mm
- Article No. 69577501



E 735/1 Mobile Unit

- With 3 levels for various inserts
- 2 built-in spray arms
- Magnetic coding strip for automatic mobile unit sensor (optional)
- Includes 3 E 750 units to secure and position inserts for optimal cleaning results
- Clearance: (Level 1 = Lower Level)
1: 203 mm H, 482 mm W, 590 mm D
2: 203 mm H, 488 mm W, 546 mm D
3: 133 mm H, 488 mm W, 546 mm D
- Dimensions: Height 522 mm
Width 530 mm
Depth 600 mm
- Article No. 69573502



E 701/1 Mobile Unit

- With 4 levels for various inserts
- 3 built-in spray arms
- Magnetic coding strip for automatic mobile unit sensor (optional)
- Includes 3 E 750 units to secure and position inserts for optimal cleaning results
- Clearance: (Level 1 = Lower Level)
1: 87 mm H, 482 mm W, 590 mm D
2/3: 87 mm H, 488 mm W, 546 mm D
4: 223 mm H, 488 mm W, 546 mm D
- Dimensions: Height 461 mm
Width 530 mm
Depth 600 mm
- Article No. 69570102



E 757 Mobile Injector Unit

- For 1 - 6 large items (ex. carboys)
- 6 injector jets with plastic star bases
- Height adjustable holding frame with 6 short and 4 long struts for various glass diameters
- Magnetic coding strip for automatic mobile unit sensor (optional)
- Injector Dimensions: Height 10 mm
Width 1.5 mm
Depth 300 mm
- Unit Dimensions: Height 346 mm
Width 530 mm
Depth 600 mm
- Article No. 69575701



E 702 Modular Unit

- For use with E 701, 1 built-in spray arm
- Creates removable 5th level
- Includes 1 E 750 unit to secure and position inserts for optimal cleaning
- Clearance with E 701:
4: 87 mm H, 488 mm W, 546 mm D
5: 81 mm H, 488 mm W, 546 mm D
- Dimensions: Height 160 mm
Width 530 mm
Depth 560 mm
- Article No. 69570201



Modular Inserts: Direct Injection



Modular Baskets: Design Your Own

Miele's modular basket concept allows the user to design a unique wash basket through the combination of various inserts, connected by a simple docking system.


When configured, the wash basket can be a combination of only injection cleaning, only non-injection cleaning, and any variation in-between.

For even greater throughput, the modular basket configuration provides easy load, unload, and transfer. Unlike other manufacturers, the G 7825 modular basket allows the operator to remove one insert at a time, eliminating the need to load and unload at the machine or lift the entire large, heavy basket.



E 741/1 Modular Mobile Unit

- Accepts 3 levels of modular inserts
- For single, double, or 3 level cleaning
- Injection, standard, or combo cleaning
- Magnetic coding strip for automatic sensing of a specific wash routine
- Two top level tracks allows flexibility per wash load
- Dimensions: Height 680 mm
Width 530 mm
Depth 600 mm
- Article No. 69574102

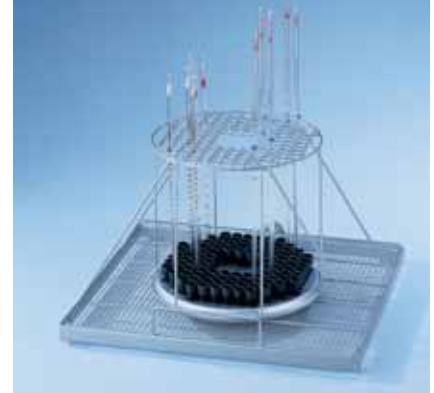

E 745/1 Modular Injector Insert

- Injector unit for 104 pipettes
- Each section is 16 mm x 16 mm
- Forced-air drying through injectors
- Frame supports wide mouth glassware
- Dimensions: Height 288 mm
Width 492 mm
Depth 496 mm
- Article No. 69574502



E 742 Mobile Unit

- Support frame with spray arm
- Standard cleaning frame for levels 2, 3, or 4
- Supports two half inserts such as the E 106 of AK 12
- Dimensions: Height 112 mm
Width 492 mm
Depth 496 mm
- Article No. 69574201




E 746 Modular Injector Insert

- Injector unit for 23 pipettes
- Holds 10 pipettes up to 560 mm long
- Holds 13 pipettes up to 490 mm long
- Forced-air drying through injectors
- Frame supports wide mouth glassware
- Allows a second cleaning level
- Dimensions: Height 330 mm
Width 492 mm
Depth 496 mm
- Article No. 69574601



E 755 Modular Injector Insert

- For narrow neck items, 25-100 ml
- Injectors removable for customization
- Forced-air drying through injectors
- 36 injectors (E 470), 2.5 x 90 mm
- Dimensions: Height 130 mm
Width 492 mm
Depth 496 mm
- Article No. 69575501

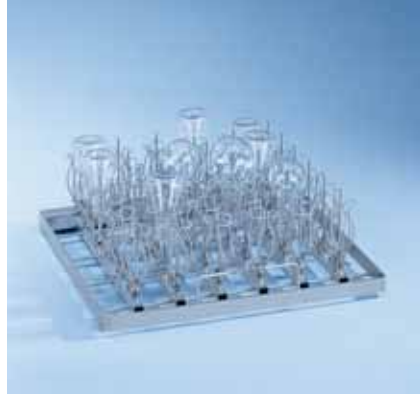

E 743 Modular Injector Insert

- For narrow neck items (up to 500 ml)
- Injectors removable for customization
- Forced-air drying through injectors
- 36 injectors (E 351), 4 x 160 mm
- 36 Mielava injector clips (E 353) for securing glassware in place
- Dimensions: Height 190 mm
Width 492 mm
Depth 496 mm
- Article No. 69574301



E 752 Modular Injector Insert

- For narrow neck items, 100 - 1000 ml
- 25 injectors (12 x E 351 jets 4 x 160 mm and 13 x E 352 jets 6.0 x 220 mm)
- 25 Mielava injector clips (12 x E 353 and 12 x E 354) for securing glassware in place
- Dimensions: Height 260 mm
Width 492 mm
Depth 496 mm
- Article No. 69575201



E 744 Modular Injector Insert

- For narrow neck items, 500-1000 ml
- Injectors removable for customization
- Forced-air drying through injectors
- 16 injectors (E 352), 6 x 220 mm
- 16 Mielava injector clips (E 354) for securing glassware in place
- Dimensions: Height 250 mm
Width 492 mm
Depth 496 mm
- Article No. 69574401



E 747 Modular Injector Insert

- For use with E 741 mobile unit
- 104 injectors, 2.5 x 110 mm
- Ideal for cleaning centrifuge tubes, vials, test tubes, or other small items
- Forced-air drying through injectors
- Frame supports wide mouth glassware
- Dimensions: Height 168 mm
Width 492 mm
Depth 496 mm
- Article No. 69574701

Baskets and Inserts: For Wide-Mouth Glassware



E 742 Mobile Unit

- Support frame with spray arm
- Standard cleaning frame for levels 2, 3, or 4
- Supports two half inserts such as the E 106 of AK 12
- Dimensions: Height 112 mm
Width 492 mm
Depth 496 mm
- Article No. 69574201



E 484 Stainless Steel Basket

- Insert 1/1, basket version
- E 486 Stainless Steel Support available for tall objects such as trays
- Dimensions: Height 65 mm
w/ handle 150 mm
Width 470 mm
Depth 480 mm
- Article No. 69548401

E 486 Stainless Steel Support

- Article No. 69548601



E 109 Half Insert

- For 21 beakers (up to 250 ml) or other glassware
- 3 rows of 140 mm tall paired holders
- For upper, lower, or half injector baskets
- Dimensions: Height 155 mm
Width 230 mm
Depth 460 mm
- Article No. 69510901

E 144 Half Insert

- For 18 beakers (up to 250 ml)
- 3 rows of 120 mm tall paired holders
- For upper, lower, or half injector baskets
- Dimensions: Height 131 mm
Width 200 mm
Depth 445 mm
- Article No. 69514401

E 110 Half Insert

- For 10 beakers (250 to 600 ml) or other glassware
- 2 rows of 160 mm tall paired holders
- For upper, lower, or half injector baskets
- Dimensions: Height 175 mm
Width 230 mm
Depth 460 mm
- Article No. 69511001

E 111 Half Insert

- For 8 beakers (600 to 1000 ml) or other glassware
- 4 rows of 190 mm holders (24 total)
- For use in lower or half insert baskets
- Dimensions: Height 205 mm
Width 230 mm
Depth 460 mm
- Article No. 69511101



E 106 Half Insert

- For lower or half injector baskets
- 26 total spring clips to secure glassware
16 small spring clips, 105 mm Tall
10 large spring clips, 175 mm Tall
- Distance between clips ~60 mm
- Dimensions: Height 186 mm
Width 195 mm
Depth 430 mm
- Article No. 69510601



E 106/1 Half Insert

- For upper, lower, or half injector baskets
- 26 small spring clips to secure glassware, 105 mm Tall
- Distance between clips ~60 mm
- Dimensions: Height 116 mm
Width 195 mm
Depth 410 mm
- Article No. 69510602




E 106/2 Half Insert

- For lower or half injector baskets
- 13 large spring clips to secure glassware, 175 mm Tall
- Distance between clips ~85 mm
- Dimensions: Height 186 mm
Width 180 mm
Depth 420 mm
- Article No. 69510603



AK 12 Half Insert Basket


- Half insert basket with handles, for beakers, bottles, and various other glassware which do not require direct injection; insert also ideal for other labware (funnels, stirrers, stopcocks)
- Dimensions: Height 67 mm
w/ handle 127 mm
Width 225 mm
Depth 442 mm
- Article No. 69501201

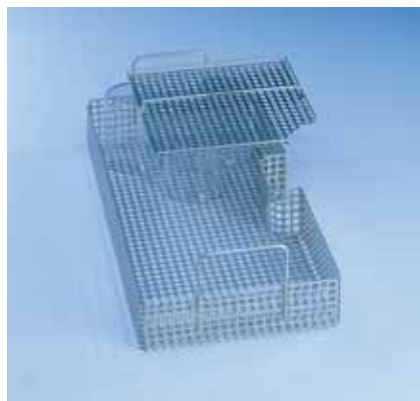

A 14-06 - 1/4 Lid (for AK 12, 2 required)


- Shown above with AK 12 Insert Basket
- Article No. 69745002

Machine Accessories

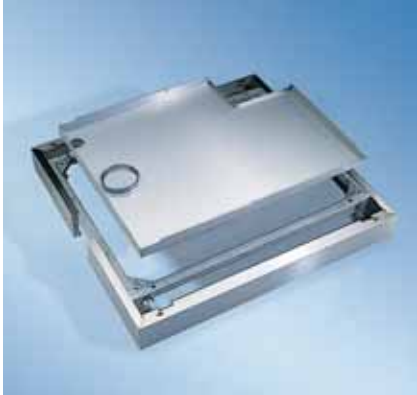

SBW/2 Base

- Stationary base
- Simplifies machine installation
- Acts as a drip tray
- Cut-out for possible floor drain
- Dimensions: Height 100 mm
Width 900 mm
Depth 734 mm
- Article No. 69371002

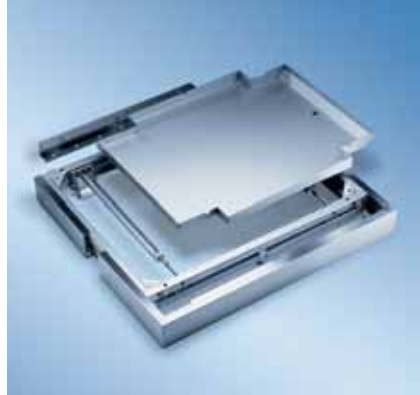

SBWR/2 Roller Base

- Mobile base equipped w/ locking rollers
- Allows machine to be easily moved away from wall for service access
- Simplifies machine installation
- Recommended if machine must be installed <18" from wall
- Acts as a drip tray
- Dimensions: Height 100 mm
Width 900 mm
Depth 734 mm
- Article No. 69371004



MF/3 Transfer Cart

- For safe and easy transfer of baskets
- Unique patented design suspends mobile unit in the air, allowing the cart to load from the front or side, saving valuable floor space
- Dimensions: Height 1182 mm
+/- 100 mm
Width 660 mm
Depth 807 mm
- Article No. 69200107



Indicator Light Box

- Indicates progress of wash cycle

3 Light Status Indicator Box

- Shown above
- Article No. PROF18975T999

5 Light Status Indicator Box

- Article No. 78270022



DK 25/26 Steam Condenser

- Requires MAV 25/26 Steam Condenser Enclosure
- Article No. 69230005

MAV 25/26 Steam Condenser Enclosure

- Required with DK 25/26 Steam Condenser
- Article No. 69210008

ML Magnetic Code Strip

- Article No. 69745702



PRT/1 Printer

- Shown above
- Article No. 69221102

External Printer

- Article No. 5035-033-01

G 7825/G 7826 Trim Kit

- For in-wall installations
- 3" wide stainless steel trim to cover wall opening most commonly for pass-through installations
- Article No. PROFTRIMKITG7825

Machine Accessories: G 7895/1 Pure Water System



G 7895/1 Pure Water System

- Stainless steel (shown) or white cabinet
- Inc. conductivity meter (in microS/cm)
- Inc. (1) E 318 cartridge & (1) E 315 resin
- Unit holds max. of two E 318 cartridges
- Dimensions: Height 850 mm
w/out lid 820 mm
Width 300 mm
Depth 600 mm
- Two pressure hoses, approx. 4' long
- Water intake to 3/4" Male Hose Thread
- Min. 36 psi incoming water pressure

G 7895/1 - Stainless Steel

- Article No. 69789504

G 7895/1 - White

- Article No. 69789503



E 318 Stainless Steel Cartridge

- Stainless steel cartridge
- Pressure-proof w/ vent & relief valve
- Holds 20 Liter of E 315 disposable resin
- Diameter 240 mm, Height 570 mm
- Article No. 69431801



E 316 Refill Kit for Resin Replacement

- Refill kit for resin replacement
- Inc. 30 Liter plastic barrel and funnel
- Capacity: 12 gallon/45 Liter
- Article No. 69431601



E 315 Disposable Resin

- Disposable resin, 20 Liter
Two 10 Liter bags
- Vacuum-packed plastic bags
- Inc. filter bag for replacement
- Article No. 69431501



Service and Validation



Validation Documentation

- Complete IQ/OQ
- User-friendly, step-by-step check-offs
- Covers all alarm conditions
- Assures safe validation
- Assures documents are up-to-date



Validation Execution

- Miele-direct technicians
- Very professional and expertly trained
- Prompt
- Set, competitive pricing alleviates cost overruns



Preventive Maintenance Plans

- Assures you of maximum machine performance and longevity
- Various plans to meet your requirements and budgets
- Performed by Miele direct or factory trained technicians

Miele
PROFESSIONAL

G 7825

PROGRAM
PROGRAMME-SUP
DOOR OPEN

I-O



ME 1:MINI
RUEY
>PROGRAMMING< [D]



PC ●

Miele Professional

9 Independence Way
Princeton, NJ 08540
USA

& 800.991.9380

☎ 609.419.4241

🌐 labwasher.com

Miele Limited

161 Four Valley Drive
Vaughan, ON L4K 4V8
Canada

& 800.643.5381

☎ 905.532.2290

🌐 miele.ca

Miele S.A. de C.V.

Av. Santa Fe 170, German Centre: 0-4-2
Col. Lomas de Santa Fe
C.p. 01210. Mexico, D.F.

& (+5255) - 8503 9870 al 73

☎ (+5255) - 8503 9874

🌐 miele-professional.com.mx

Every effort has been made to ensure that the information given in this brochure is correct at the time of printing. However, due to continuous product improvement, Miele reserves the right to make changes to products and technical data without prior notice.